



PathoSense

2024

CONFIDENTIAL

The content of this document is confidential and intended for the training of employees of official PathoSense Partnerlabs. It is strictly forbidden to share any part of this document with any third party, without a written consent of PathoSense.



Wet lab flow diagnostics PathoSense

WET LAB VI - RUN SET-UP

**Set-up of a sequencing run on the MinKNOW software
and generation of a sequencing run file**

Timing ~ 5 min

1. Open the MinKNOW software and click on **'Start sequencing'**
2. Fill in the **experiment name**
3. Click on the position of the correct flow cell
4. Copy the **experiment name** and paste it in the Sample ID
5. Click on **'Continue to kit selection'**
6. Select the used kit (Rapid Barcoding Kit 24, **SQK-RBK114.24**)
7. Click on **'Continue to run options'**
8. Click on **'Options'**, change **'Run time'** to **'9' Hrs** and click on **'Apply rules'**
9. Click open **'Advanced options'**, turn **off 'Reserve pores'** and change **'Time between pore scans'** to **'1' Hours**

10. Click on '**Continue to analysis**'
11. In the section 'Basecalling', click on '**Edit options**', change '**Basecalling model**' to '**Super-accurate basecalling**' and click on '**Apply**'
12. In the section 'Barcoding', click on '**Edit options**', turn on '**Trim barcodes**' and click on '**Apply**'
13. Click on '**Continue to output**'
14. Select the correct output location (**PathoSense_Diagnostics**)
15. Click on '**Options**', change '**Qscore**' to '**1**' and click on '**Apply**'
16. Click on '**Continue to final review**'
17. Click on '**Start**'
18. While the flow cell is getting to temperature, go to cloud.pathosense.com and log in with your credentials.
19. On the Runs page, make a new run by setting the "**Size of run**" to the correct size and pressing "**+ New run**"
20. Select the run by clicking anywhere on the run and add the samples by clicking them
***Note:** only arrived samples can be added, if the samples are in "En route", set the analysis to "Arrived" on the analysis page*
21. Click the green **save icon** and select the used ONT barcode for each of the samples
22. Press "**Download**" to download the run file
23. Press the blue "**Start run**" button to move the run to "**Running**"
24. When the flowcell has reached it's temperature, the sequencing output folder should appear in the "**PathoSense_Diagnostics**" folder
25. Copy the downloaded run file and paste it in the sequencing folder, to the same directory that contains the "**fastq_pass**" and "**pod5_pass**" folders

